



Improved digestions for better analysis



BD50 & BD28 SPECIFICATIONS

POWER REQUIREMENTS

Block: 220-260VAC, 50/60 Hz
Controller: 100-260VAC, 50/60 Hz

OPERATION REQUIREMENTS

Digestion block must be operated in a fume hood. Controller should be operated outside of fume hood.

FOOTPRINT

Block: 60 cm (23.5") W x 28 cm (11") D
Controller: 25 cm (10") W x 24 cm (9.5") D

TEMPERATURE RANGE

Ambient to 450°C

SET POINT ACCURACY

+/- 2°C

CAPACITY

50 x 75ml tubes, 50 x 100ml tubes, 28 x 250ml tubes

SYSTEM CONFIGURATIONS

DG-U-A021 BD28s system (28 place) consists of a Heating Block, Programmable Controller and Tube rack/draftshield

DG-U-A024 BD50s system (50 place) consists of a Heating Block, Programmable Controller and Tube rack/draftshield

ACCESSORIES

- 4003035** Digestion tubes, glass, 75ml, straight (50pk)
- 4003005** Digestion tubes, glass, 75ml, volumetric (50pk)
- 4003036** Digestion tubes, glass, 100ml, straight (50pk)
- 4003006** Digestion tubes, glass, 100ml, volumetric (50pk)
- 4003031** Digestion tubes, glass, 250ml, straight (28pk)
- 4003001** Digestion tubes, glass, 250ml, volumetric (28pk)
- 5001241** Cooling stand for 28 place tube racks
- 5001511** Cooling stand for 50 place tube racks
- 8027245** Teardrop stopper, glass (75/100ml tubes) for enhanced reflux
- 8029292** Teardrop stopper, glass (250ml tubes) for enhanced reflux
- 8027565** Splash-head, glass, for 250ml tubes
- 5001231** Tube rack/draftshield, 28 place for 250ml tubes
- 8028114** Tube rack/draftshield, 50 place for 75ml or 100ml tubes



www.seal-analytical.com

SEAL Analytical is a global company with offices worldwide - contact us at:

SEAL Analytical, Inc.

Mequon Technology Center
10520-C Baehr Rd.
Mequon, WI 53092
United States
Tel: +1 (262) 241 7900
Fax: +1 (262) 241 7970
sales@seal-us.com

SEAL Analytical Ltd.

2 Concorde Close
Segensworth
Fareham
Hampshire PO15 5RT
United Kingdom
Tel: +44 (0) 1489 864400
Fax: +44 (0) 1489 880531
mail@seal-analytical.com

SEAL Analytical GmbH

Werkstrasse 5
D-22844 Norderstedt
Germany
Tel: +49 (0)40 60 9292 9-00
Fax: +49 (0)40 60 9292 9-02
info.germany@seal-analytical.com

SEAL Analytical Shanghai

Fortune International Plaza
Silver Building, Room 614
No. 43, Guo Quan Road
Shanghai 200433
China
Tel: +86 21 3362 5002
Fax: +86 21 3362 5002



BD50 & BD28

BLOCK DIGESTION SYSTEM


BD50 & BD28




28 place block with digestion tubes in the heating position



50 place block with digestion tubes in the cooling position

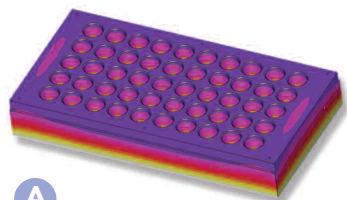

A
The BD50s and BD28s unique flat plate heating design provides even heating leading to better quality digestions.

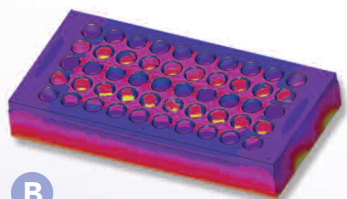

B
Common rod heating blocks have large temperature variations across the block resulting in poor, uneven digestions.

BETTER PRODUCTIVITY

The BD50s and BD28s Programmable Digestion Systems perform acid digestion of large numbers of samples under controlled conditions and are available in two different formats:

- ▶ 28 place for 250ml digestion tubes
- ▶ 50 place for both 75ml and 100ml digestion tubes

Designed for durability and robustness, a BD50s/BD28s digestion system includes a block digestion unit, microprocessor controller, tube rack/draftshield, a set of digestion vessels and a cooling stand that supports the tube rack above the block.

If you are looking to automate your digestions, the BD50s and BD28s offers enhanced productivity, better quality digestions and increased safety.

MORE RELIABLE DIGESTIONS

The BD50s and BD28s Digestion Systems heat samples at a controlled rate, in the presence of one or more acids, to the required temperature then holds the samples at that temperature until digestion is complete.

The BD50s and BD28s' unique flat plate design results in an even temperature distribution across the block and more precise temperature control. The temperature of each digestion position in the BD50s and BD28s is tightly regulated ensuring all samples in the batch are digested equally (diagram A).

Other block designs use a more common rod heater design that results in a greater temperature variation across the digestion block (diagram B).

The different coloured zones shown on the BD50s and BD28s block illustrate the temperature isotherms across the block at 400°C.

In addition, your digestion methods are stored electronically to provide reproducibility from batch- to-batch, operator-to-operator. You can store up to 20 programs, each with up to 30 steps, allowing you to automate the most complex digestion procedure.

- ▶ **As a market leader, SEAL has over 1,000 applications available and under continual development. Markets include water, wastewater, soil, plant, fertilizer, food and beverage. Please contact us for your specific application.**

SAFETY

The excellent temperature stability of the BD50s and BD28s eliminates the need to move individual samples to a more "favourable position" during the digestion.

Controlled heating rates, as low as 1°C per minute, protect against bumping and splattering of reactive samples.

The BD50s and BD28s Controller senses power failures and will record the digest conditions and hold the digest to protect your samples. An in-built safety cut-out protects from overheating. The Controller is separate from the Block protecting the electronics and the operator from acid fumes.

IDEAL FOR

- ▶ Total Kjeldahl Nitrogen
- ▶ Total Phosphorus
- ▶ Metals
- ▶ Mercury
- ▶ C.O.D.
- ▶ and more digestions

ECONOMICAL

Apart from power and reagents, the BD50s and BD28s have no ongoing consumable costs and do not use any components with a limited service life.


The provision to program an elevated standby temperature for rapid start up enables you to process more batches per day. In addition, you can safely run a digest overnight so you have a full batch of digested samples ready for analysis at the beginning of each day. Automatic operation frees key laboratory personnel from the tedium of managing long digestion processes and provides an early start for analysis.

The BD50s and BD28s also have a small footprint to save you valuable fume cupboard space.

ACCESSORIES



Tube rack/draftshield
Enables all tubes to be easily transported and keeps the radiated heat around the tubes during digestion.

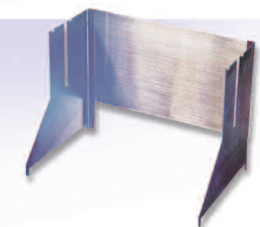

Cooling Stand
Supports the tube rack above the block to facilitate rapid cooling.



Digestion Tubes
Available in straight sided tubes or as volumetric tubes with a constricted neck and class "A" calibration mark.



Teardrop stoppers
Captures and recondenses digest fumes reducing acid loss and preventing cross contamination.



Splash-head
For safe reagent addition during a digestion to 250mL digestion tubes.



PLEASE VISIT

www.seal-analytical.com

or contact your local SEAL Analytical representative about your specific application.